

## 5.0.5244

- *New Feature* – User based Time Zone and Date Settings
- *New Feature* – Quick Group Selection on Reports
- *New Feature* – 4hr Sensor Summary Report
- *New Feature* – Current Readings Report
- *New Feature* – VFC Report
- *New Feature* – Global Acknowledgement
- *New Feature* – Calibration Documentation
- User Security Improvements
- Reports Enhanced
- UI Updated
- Active Directory Integration Enhanced
- Collector Service Updated
- Database Updated
- Installation Package Modified
- Popup Utility Updated
- EZLink Updated

### 5.0.4316

- **New Feature** – Support for new Digital Steam trap and Dual Input Sensors
- Collector Service Updated
- Notification Service Updated
- Installation Package Modified
- Sensor Registration Improved
- Reports Enhanced

### **5.0.4170**

- Collector Service Updated
- Updated SSRS Reports
- Tags (Search Filter) Updated

### 5.0.4139

- **New Feature** – Windows 8 / 2012 + SQL 2012 Support
- **New Feature** – Reports now have selectable date ranges
- **New Feature** – Acknowledge / Clear from Sensor Details Page
- **New Feature** – Acknowledge Alarms Only Permission
- **New Feature** – Extended Data Retention Options
- **New Feature** – Sensor Groups can be alphabetized
- Improved Sensor Calibrations
- Collector Service Updated
- Updated SSRS Reports
- User Interface Updated
- SQL Express Capacity Warning Improved
- Change Sensor Types Updated
- TempTrak Login Improved

## 5.0.3338

- **New Feature** – TempTrak now Cluster-Aware
- **New Feature** – Cart Reports
- **New Feature** – SMS Groups (International Customers)
- **New Feature** – Tags (Search Filter)
- **New Feature** – Extended Logging Profiles (12hr/24hr)
- Improved Sensor Calibrations
- Collector Service Updated
- Notification Service Updated
- Enhanced detailed sensor information
- Sensor Topology updated
- Updated SSRS Reports
- Improved Map Layout icons

### **5.0.3255**

- Updated EZLink
- Improved support for long Group and Sensor names
- Updated built-in help content
- Enhanced view creation
- Updated SSRS Reports
- Current Alerts updated with Grouping and Sensor Description
- Sensor Attributes updated with Grouping and Sensor Description
- Fixed issue with change sensor type transmission interval
- Collector Service Updated

### 5.0.3186

- **New Feature** – International Report Compatibility
- **New Feature** – Sensor History & 24hr History Chart now show average reading
- **New Feature** – Internet Explorer 10 Support
- **New Feature** – NIST Validation now supported on Dual Input 900MHz Transmitters
- Enhanced profile deletion display prompt
- Improved sensor delete message prompt
- Fixed dropdown display issue with notification groups
- Resolved sensor name display issue in EZLink
- Improved multiple database login page
- NIST Validation Report updated to display all validated probes
- Updated sensor reading charts
- Enhanced Global Search
- Updated CSS for better browser support
- Collector services updated
- Updated SSRS Reports

**5.0.3141**

- Database Update script modified for greater SQL compatibility



## 5.0.3126

- **New Feature** – Sensor Alarm limits are now displayed on SSRS Sensor History Report
- **New Feature** – SQL Express Database Size limit warning now displayed on login
- **New Feature** – User Access Permissions have been separated out for specific profile types
- **New Feature** – Enabled support for direct web interface access on multiple database installations
- Fixed display issue with repeaters on current readings
- Collector Service updated for Legacy DocuTemp transmitters
- Updated EZLink to resolve issue where 24hr History would not display for all sensors
- Changed Report Access Permissions to allow users to select the run option on SSRS Reports
- Improved Group Audit Charts allowing anyone with Reports access the ability to run the Report
- Resolved issue with Celsius conversion on Group Audit Chart SSRS Report
- Corrected issue when registering Wi-Fi Dual Temperature transmitters
- Improved Update Files script for software upgrades
- Enhanced IE7 support
- Fixed display issue with deleted notification profiles
- Updated security roles
- Fixed logging profile enabled flag display issue
- Enhanced notification service support for dynamic notification types

### 5.0.3085

- **New Feature** – Support for new Wi-Fi 802.11 b/g/n Transmitters
- Removed duplicate probe ID from alarm notification data
- Corrected issue with Daylight Savings Time in SSRS reports and User Interface
- Updated sensor audits in view details to indicate if no audit performed
- Improved collector restart warning message when performing sensor replace and relocate
- Enhanced transmitter registration ID Validation
- Adjusted timeout on sensor delete function
- Fixed report scheduler issues with midnight report times
- Resolved issue where alarm profiles were not correctly converting to Celsius
- Updated notification groups to allow blank groups
- Changed time validation in profile configuration
- Fixed issue with sorting on Sensor ID
- Corrected issue with blank custom message formats
- Updated profile associations so they are removed upon deletion
- Updated User Settings to display selected options upon users first login
- Resolved permissions issue with sensor history chart and NIST SSRS reports
- Removed extra date information from Alerts by Day report

### **5.0.3042**

- Updated menu cache for custom report list to improve performance
- Resolved problem with transmitter replacement case sensitivity
- Changed reporting periods for Daily Summary Report so it would function with the users selection
- Corrected issue with default reporting period in user settings
- Fixed issue that caused group duplications in details view
- Updated Topology Report to highlight repeaters
- Updated Group/View Map sensor icons to show if sensor is in alarm range
- Removed extra probe selection option when registering DocuTemp Analog Transmitter
- Updated help descriptions
- Changed scale on contact transmitter charts
- Improved loading of sensor details
- Corrected issue with SQL 2005 Reports preventing reports from displaying

## 5.0 RTM

- **New Feature** – Dashboard view: Provides high level overview
- **New Feature** – EZLink: Providing users with a mobile interface for clearing alarms
- **New Feature** – Scheduled reporting (not available when using SQL Express)
- **New Feature** - Custom sensor ordering: Users can order sensors in groups / groups within view
- **New Feature** – Sorting / Filtering / Searching / Paging features for greater usability
- **New Feature** – Help documentation is specific to the current page the user is viewing
- **New Feature** – Ability to add Custom Reporting
- **New Feature** - Reports generated using SSRS (SQL Server Reporting Services)
- Updated user interface
- Removed Java requirement from web interface: java is not needed to run TempTrak
- Removed ActiveX Components from web interface
- Removed Auto-Registration of hardware
- Enhanced user security allowing for more granular security rights
- Enhanced profile security allows for easily assigning profiles to specific views
- Enhanced sensor security registration
- Enhanced sensor reading charts